

Development of Trace Carbon Dioxide Isotopologue Analyzer - Performance Evaluation Study

BJ Clark¹, Graham Leggett¹, Doug Lynch¹, Mark Johnson¹, Anatoly Komissarov¹, Israel Begashaw¹

1. LI-COR Biosciences, Lincoln, Nebraska, USA.

Introduction

We report on the development and performance evaluation of a soon-to-be released trace gas analyzer for the measurement of atmospheric $\delta^{13}\text{C}$, $\delta^{18}\text{O}$, $\delta^{17}\text{O}$, ^{13}C , ^{12}C , ^{18}O and ^{17}O of CO_2 . Carbon Dioxide isotopes can help answer many questions, such as emission source attribution and, via their use as a tracer, to examination of naturally occurring processes in plants or soil. Based on Optical Feedback Cavity Enhanced Absorption Spectroscopy (OF-CEAS), the analyzer offers both the sensitivity and stability necessary for long-term atmospheric monitoring, DIC ocean measurements, soil, plant physiology and mud gas isotope logging applications.

Analyzer Specifications



- Total Weight: 10.5 kg (including batteries)
- Battery Life: 8 hours typical with 2 batteries
- Operating temperature range -25 °C to +45 °C

LI-7825 CO₂ Isotope Analyzer

CO₂ Measurements:

Measurements: Range: 50 to 2,000 ppm
Precision (1 σ): 0.027 ppm with 5-minute averaging
Maximum drift: <0.4 ppm per 24-hour period

$\delta^{13}\text{C}$ Measurements:

Precision (1 σ): 0.02 ‰ with 5-minute averaging
Maximum drift: <0.5 ‰ per 24-hr period

$\delta^{18}\text{O}$ Measurements:

Precision (1 σ): 0.06 ‰ with 5-minute averaging
Maximum drift: <2 ‰ per 24-hr period


$\delta^{17}\text{O}$ Measurements:

Precision (1 σ): 0.2 ‰ with 5-minute averaging
Maximum drift: <6 ‰ per 24-hr period


Test Setup

Data for the Allan Deviation plots were collected over a 10-day period, where, prior to the measurement of test gas, the LI-7825 prototype was powered on/warmed up while sampling ambient air for 24 hours. It was then connected to a 400 ppm CO_2 tank via stainless steel tubing for the test.


$\delta^{13}\text{C}$ Allan Deviation

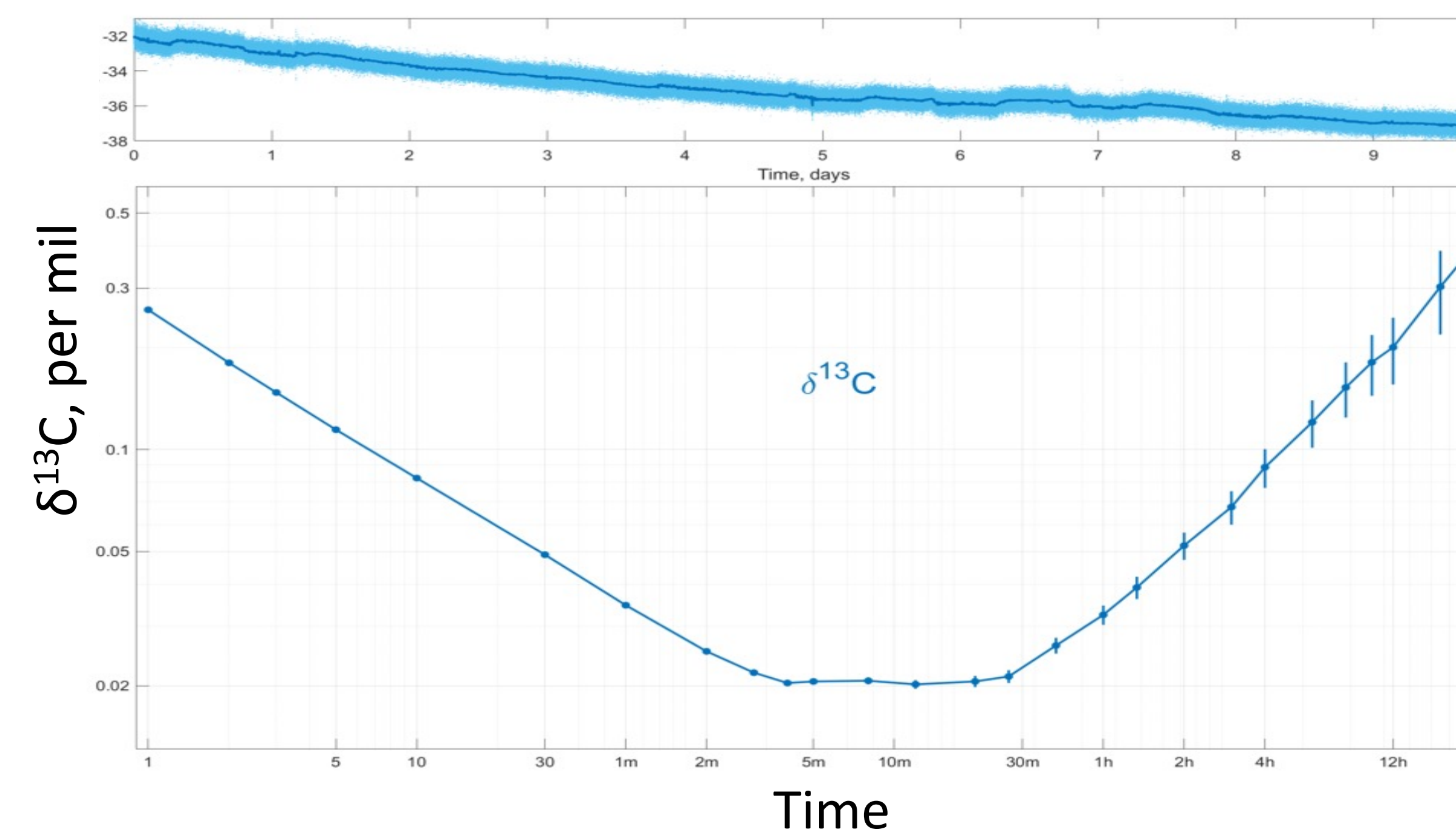


$\delta^{18}\text{O}$ Allan Deviation

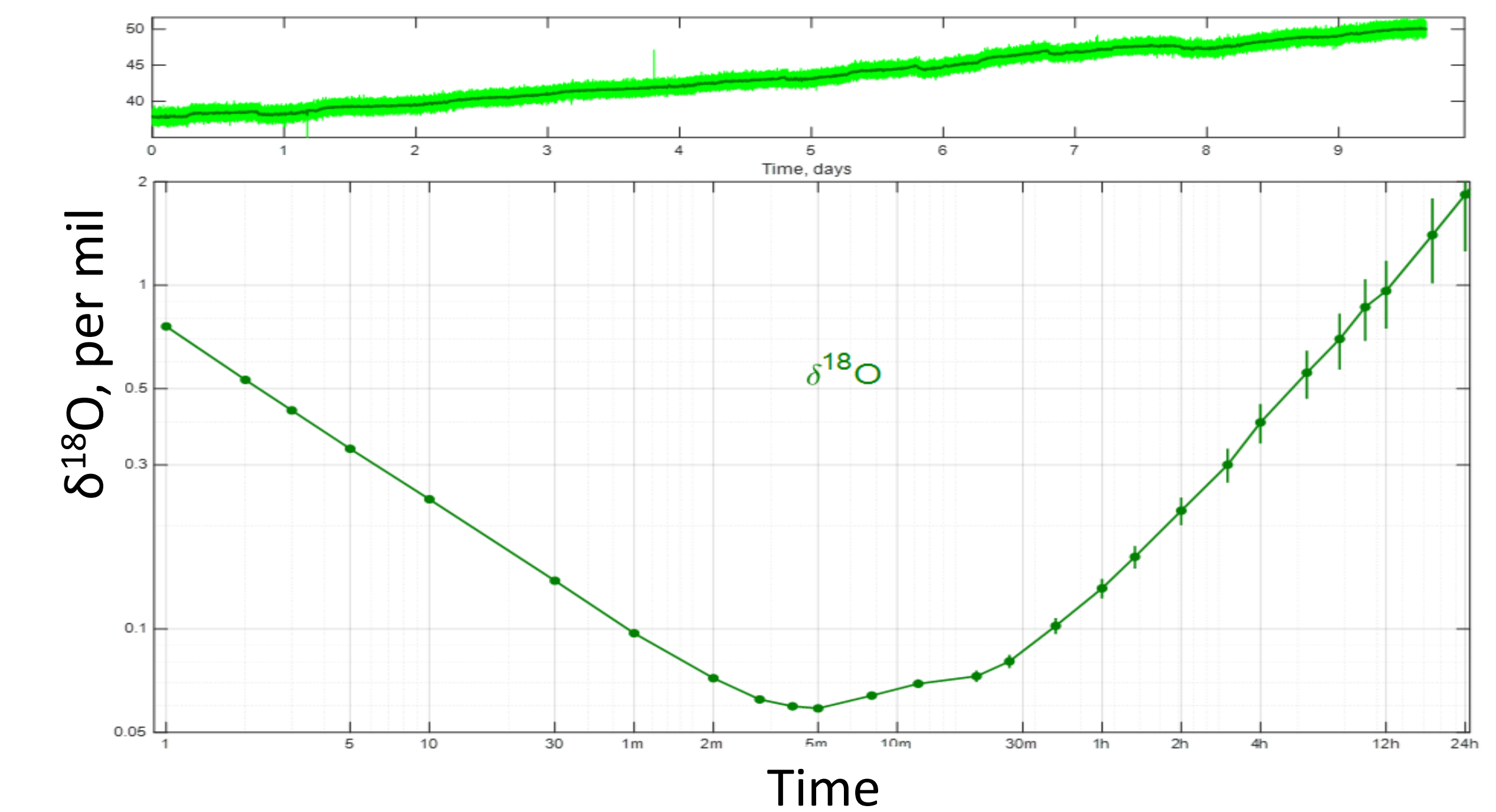

$\delta^{17}\text{O}$ Allan Deviation

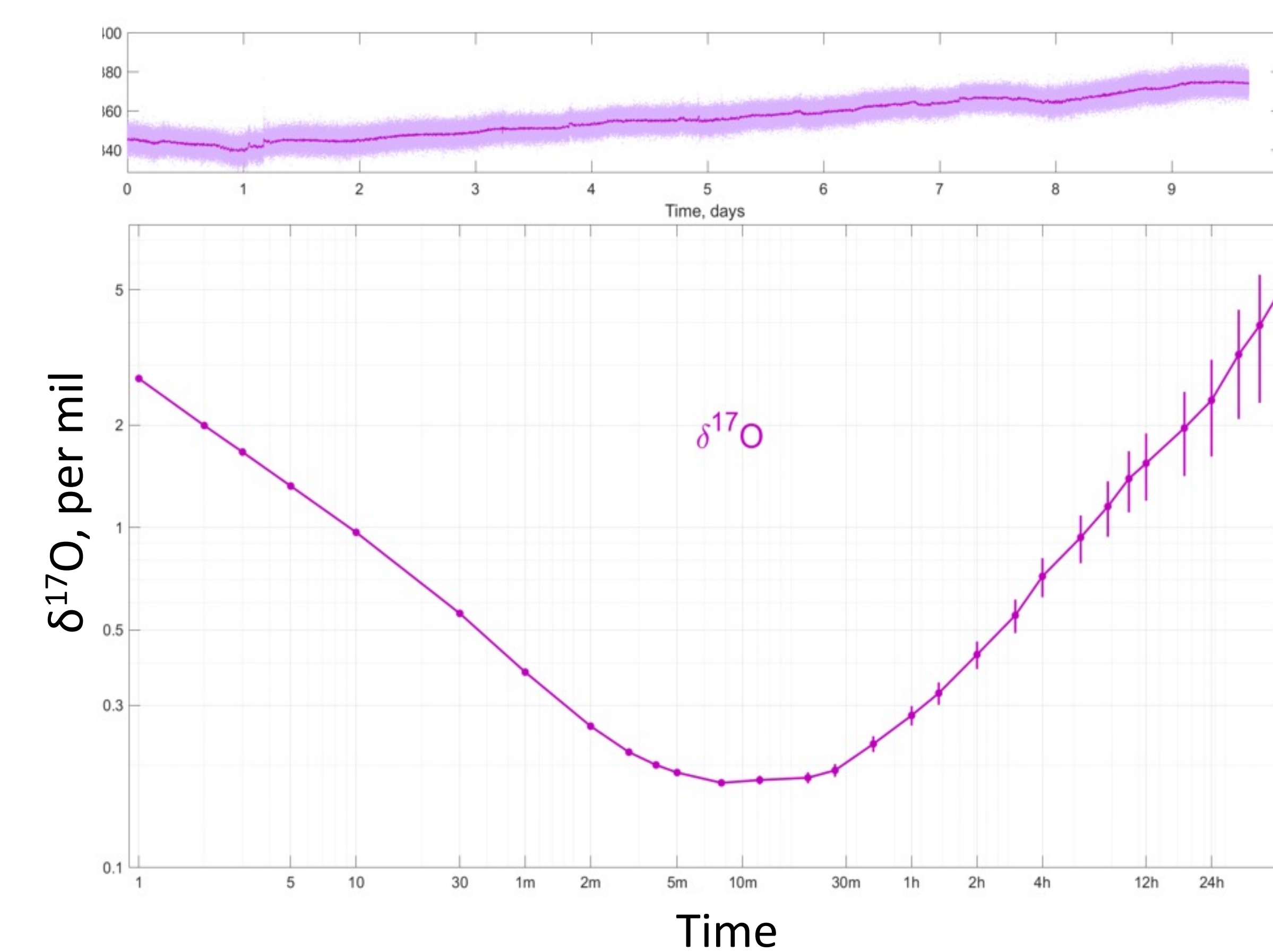

CO₂ Total Allan Deviation

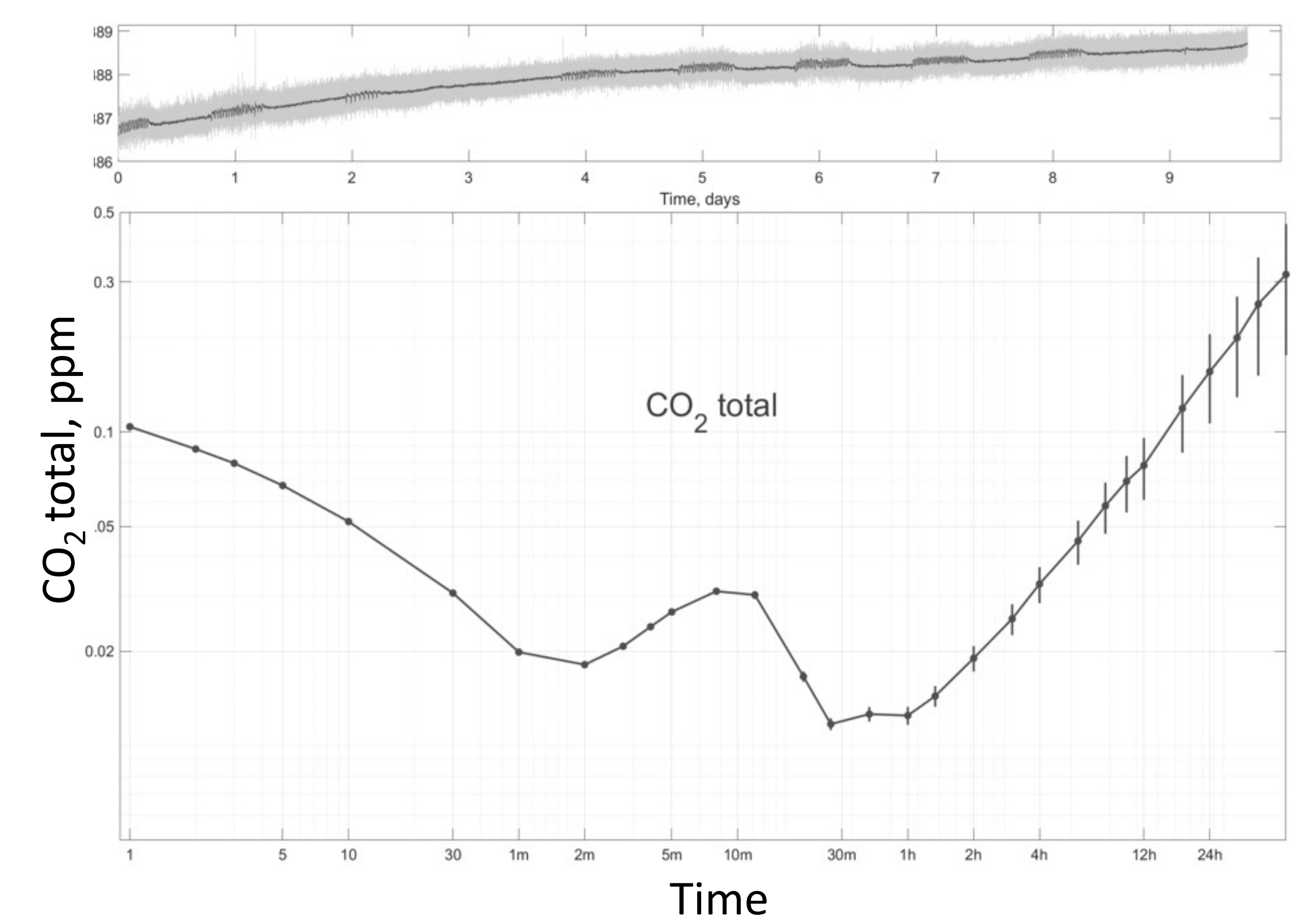

Conclusions

In conclusion, this portable and rugged instrument meets or exceeds requirements for both long-term atmospheric background measurements and offers a versatile platform for a range of mobile and agile measurements relevant to the better understanding of greenhouse gas emissions from anthropogenic and natural sources.

Acknowledgements

The initial principles and some portions of the new technology presented herein were developed in part based on the grant from the MONITOR Program by the U.S. Department of Energy Advanced Research Projects Agency - Energy (ARPA-E), under award number DE-AR0000537.